Nano Gel Extraction Kit







ADVANCED MICRODEVICES PVT. LTD.

21, Industrial Area, Ambala Cantt - 133006 (INDIA) Tel: 0171-2699290, 2699471, 2699274, Fax: +91-171-2699221, 2699008 Email: info@mdimembrane.com, Website: www.mdimembrane.com

Table of Contents

| 1. | Introduction | 3 |
|-----|---|-------|
| 2. | Applications | 3 |
| 3. | Storage Conditions | 3 |
| 4. | Quality Assurance | 4 |
| 5. | Safety Information | 4 |
| 6. | Lot Release Criteria | 4 |
| 7. | Technical Support | 4 |
| 8. | Kit contents | 5 |
| 9. | Specifications | 5 |
| 10. | Principle | 6 |
| 11. | Important Points to be Considered | 6 |
| 12. | mdi Nano Gel Extraction Procedure | 7 |
| 13. | Protocol | 8-9 |
| 14. | Trouble Shooting Guide | 10-11 |
| 15. | Product Use Limitations | 11 |
| 16. | Product Warranty and Satisfaction Guarantee | 11 |
| 17. | Ordering Information | 12 |
| | | |

1. Introduction

mdi Nano Gel Extraction Kit is a quick, convenient and economical tool to obtain highly concentrated DNA fragments (70bp-4kb) after gel analysis from standard or low melt Agarose gel. The buffer system provided in the kit facilitates efficient binding of DNA fragments from solublised gel slice onto the spin column.

Washing is done with help of wash buffers in order to remove agarose gel, salts, and other impurities from DNA samples. Purified DNA is eluted in small volumes (5µl) of low salt buffer or water for variety of downstream applications.

2. Applications

- 1. Automated Fluorescent Sequencing
- 2. Radioactive Sequencing
- 3. Restriction Digestion
- 4. Ligation
- 5. Transformation
- 6. Cloning
- 7. PCR

3. Storage Conditions

As soon as you receive the kit, open it and store mdi Nano Gel Extraction spin columns at 2-8°C.

The remaining kit components can be stored at room temperature. The kit is stable for 1 year at given storage conditions. If precipitates appear in buffers, re-dissolve at 37°C for few minutes.

4. Quality Assurance

The mdi Nano Gel Extraction kit is designed for various pre-determined specifications and user requirements such as yield, purity, ruggedness, shelf life and functional convenience.

These are produced through a well defined quality management system certified by Underwriters Laboratories, USA for ISO 9001: 2008 which ensures intra lot as well as lot to lot consistency.

5. Safety Information

The buffers and the reagents may contain irritants, so wear lab coat, disposable gloves and protective goggles while working with the Nano Gel Extraction Kit.

6. Lot Release Criteria

Each lot of Nano Gel Extraction Kit is tested against predetermined specifications to ensure consistent product quality.

7. Technical Support

At mdi, customers are our priority. We will share our experiences to assist you to overcome problems in general product usage as well as customize products for special applications. We will

- * Stimulate problems, and suggest alternative methods to solve them.
- * Make changes/ improvements in our existing products/protocols.
- * Develop special new products and systems especially to satisfy your needs.

We welcome your feedback to improve our products.

8. Kit Contents

| Contents | | Quantity | Storage Temperature | | |
|------------------------|-------|----------|------------------------|-------|--|
| Spin Columns | 50 | 250 | 1000 | 2-8°C | |
| Buffer NMG | 100ml | 500ml | 2000ml | RT | |
| Buffer GW | 60ml | 300ml | 1200ml | RT | |
| Buffer GE | 15ml | 75ml | 300ml | RT | |
| Collection Tubes | 50 | 250 | 1000 | RT | |
| Hand Book | 1 | - | - | - | |
| Certificate of Quality | 1 | - | - | - | |

9. Specifications

| Capacity of column reservoir | 800µl |
|---------------------------------------|-------|
| Binding capacity of membrane (ds DNA) | 5µg |
| Recovery of DNA (70bp-4kb) | 80% |
| Maximum weight of gel slice | 400mg |
| Minimum elution volume for conc. DNA | 5µl |
| Total eluate volume | 4µl |

10. Principle

The mdi Nano Gel Extraction kit allows isolation of high ultra pure DNA fragments from solublised agarose gel which involves:

1. Capturing of DNA from solublised gel slice on spin column

In order to facilitate adsorption of DNA fragment on to the spin column, suitable pH and salt concentration are required which is achieved by addition of binding buffer 'NMG'.

2. Washing

Subsequent to DNA binding from solubilised gel slice on spin column unwanted components like dyes, salts and agarose are washed away. Washing is done by buffer 'GW'.

3. Elution

Salt concentration and pH of elution buffer is important for elution efficiency, elution occurs at basic conditions and low salt concentration in small volumes. Elution is done with buffer 'GE'.

11. Important Points to be Considered

Centrifugation

* All centrifugation steps should be carried out at room temperature at \geq 10,000 rpm in a conventional table top microcentrifuge. In case of choking of spin column, increase centrifugation time.

Washing

- * To remove residual wash buffer spin the column for 1-2 minutes at \geq 10,000 rpm.
- * Flow through should be properly discarded before additional centrifugation step otherwise it will not be removed & may inhibit enzymatic reactions.


Elution

* Elution buffer should be dispensed on to the center of column for maximum elution efficiency. If eluate volume is lower, then centrifuge for one minute extra at \geq 10,000 rpm.

Yield

DNA yield depends on volume of elution buffer and the amount of starting material which should be as per the protocol to maximize yields.

12. mdi Nano Gel Extraction Procedure

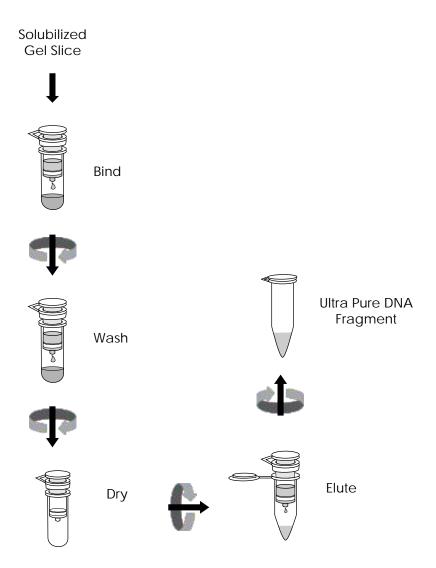

13. Protocol

- 1. Excise the DNA fragment from the agarose gel with a clean, sharp scalpel or blade. (Minimize the size of the gel slice by removing extra agarose).
- Weigh the gel slice in a colorless tube. Add 3 volumes of buffer NMG to 1 volume of gel. (100mg ~ 100µl) For example, add 300µl of buffer NMG to each 100mg of gel.
- 3. Incubate at 50 °C in a water bath for 10 minutes (or until the gel slice has completely dissolved). To help dissolve the gel, mix by vortexing the tube every 2-3 minutes, during incubation.
- After the gel slice has dissolved completely, add 1 gel volume of Isopropanol to the sample and mix by inverting the tube. For example, if the agarose gel weight is 100 mg, add 100µl isopropanol.
- Place the spin column in the collection tube in a suitable rack. To bind DNA, load the sample to the spin column and centrifuge for 1 minute at ≥10,000 rpm.

The maximum volume of the column reservoir is 800µl. For sample volumes of more than 800µl, simply load, balance micro centrifuge and spin again.

- 6. Discard the flowthrough, and using the same collection tube, wash the spin column with 500 μ l of buffer NMG by centrifuging for 1 minute at \geq 10,000 rpm.
- 7. Discard the flowthrough, and using the same collection tube, wash the spin column with 750 μ l of buffer GW by centrifuging for 1 minute at \geq 10,000 rpm.

8. Discard the flowthrough, place the spin column in the same collection tube and centrifuge for 1-2 minutes at \geq 10,000 rpm.

Important: This step is necessary to remove residual wash buffer. Residual wash buffer will not be completely removed unless the flow through is discarded before this additional centrifugation. Residual wash buffer may inhibit subsequent enzymatic reactions.

 Place the spin column in a fresh 1.5ml microcentrifuge tube (Not provided). Elute the DNA by dispensing 5µl of elution buffer GE directly on to the center of the column membrane, allow it to stand for 1 minute and centrifuge for 1 minute at ≥10,000 rpm.

Note: The elution buffer should be dispensed on to the center of the column membrane for maximum recovery. The average eluate volume is 4μ I from 5μ I elution buffer. If the eluate volume is lower, then centrifuge for one minute extra at \geq 10,000 rpm.

| 14. Trouble Shooting Guide | | | |
|---|--|--|--|
| Low or no recovery | | | |
| 1. Gel slice not solubilized | minutes. Preferably incubation should be done in a water bath instead of an oven. Use multiple columns for gel slices >400mg. Check the pH of electrophoresis buffer before use. In case it is > 8, | | |
| | Increase incubation time to 15 minutes. | | |
| | Preferably incubation should be done in a water bath instead of an oven. | | |
| 2. Gel slice >400mg | | | |
| pH of electrophoresis buffer too high | Check the pH of electrophoresis buffer before use. In case it is > 8, prepare fresh buffer. | | |
| | The electrophoresis buffer, if incorrectly prepared or used repeatedly, results in sample pH that exceeds the buffering capacity of buffer NMG and leads to inefficient DNA binding. | | |
| 4. Improper dispensing of elution buffer | Elution buffer must be dispensed on to the center of the column membrane for maximum elution efficiency. Increase incubation time 2-3 minutes more. | | |

DNA does not perform well

| Eluate contaminated with agarose | The gel slice is not properly solubilized or is more than 400mg in weight. Vortex every 2-3 minutes during incubation time. | | | | | |
|--|--|--|--|--|--|--|
| | Increase total incubation time. | | | | | |
| 2. Elution contains residual wash buffer | Spin the column for 1-2 minutes extra at \geq 10,000 rpm to remove the residual wash buffer completely. | | | | | |

15. Product Use Limitations

mdi kits are developed and manufactured for research purpose only. The products are not recommended to be used for human, diagnostics or drug purposes for which these should be cleared by the concerned regulatory bodies in the country of use.

16. Product Warranty and Satisfaction Guarantee

All mdi products are guaranteed and are backed by our

- a. Technical expertise and experience of over 30 years.
- b. Special mdi process for consistency and repeatability.
- c. Strict quality control and quality assurance regimen.
- d. Certificate of analysis accompanied with each product.

mdi provides an unconditional guarantee to replace the kit if it does not perform for any reasons other than misuse. However, the user needs to validate the performance of the kit for its specific use.

17. Ordering Information

To order please specify as below:

| Туре | | | XX | ХХ | | XX | | x | Pack Size | | |
|-----------|------|--|------|----|---|----|--|----|-----------|---------|------|
| Туре | Code | | | | | | | | Pa | ck Size | Code |
| NGEK NGEK | | | | | | | | | | 50 | 0050 |
| | | | | | | | | | | 250 | 0250 |
| | | | | | | | | | | 1000 | 1000 |
| | | | | | | | | | | | |
| Example: | | | NGEK | XX | < | ХХ | | ХХ | | Х | 0250 |

